

## **Yes, for both Windows XP and Windows 7 64 bit versions.**

First download the latest Version ProgRes CapturePro 2.7.7 from our website [www.progres-camera.com](http://www.progres-camera.com). For Login you can use for both Username and Password "Test@progres.com" ).

### ProgRes USB Camera:

- Unplug USB Camera, Install CapturePro 2.7.7, replug the USB Camera and automatically it will install the correct drivers.
- Do not forget to unplug and replug

### ProgRes FireWire Cameras

- Unplug Firewire Camera.
- Install CapturePro 2.7.7.
- Go to your CapturePro Install directory /program files / jenoptik driver and install the Thesycon T1394-Bus-Driver.

For both systems

## **Are 64 bit compatible drivers and software available for ProgRes?**

Friday, 09 April 2010 21:31

---

- Replug the camera and start the ProgRes CapturePro Software.
- Sometimes you need to restart the computer.